

About

BioClean Sterile Pharma Covers™ provide protection from environmental impurities during storage, use and transportation. They also minimize the need for autoclave tape during wrapping and save time when prepping glassware, tubing etc for autoclaving. Easy removal minimizes particulate generation and facilitates aseptic technique. Available in a range of diameters. Bespoke sizing available on request.

Specifications

MATERIAL:

Tyvek® material, IPA resistant coloured thread and high quality elastic

COMPATIBILITY:

ISO Class 4

COLOUR:

White

STERILIZATION:

Gamma irradiation

Features

- Breathable microbial barrier
- Rubber elastic provides secure seal
- Lint-free, moisture and puncture resistant
- IPA resistant coloured thread
- Form fitted
- Fast drying time
- Easy to apply and remove

LINT-FREE

MOISTURE AND PUNCTURE RESISTANT

PROVIDES SECURE SEAL



Billy Bioclean©

Quality Standards

- Processed in an NEBB certified ISO Class 4 environment
- Manufactured in a facility holding ISO 9001: 2008 certification

Sterilization

Method: Gamma irradiation
Minimum Dose: 25kGy
Sterility Assurance Level: 10^{-6}

Shelf Life & Storage

Three (3) years from date of manufacture. Store in a dry, cool place (<40°C) away from direct sunlight and fluorescent light.

Test Results

Typical Helmke Drum Test Results			
size(μm)	Average counts per minute		
	BPC-008	BPC-041	BPC-061
≥0.3	104	184	223
≥0.5	93	155	208
≥5.0	36	49	66

Endotoxin Test Results	
Method, USP <85> (Kinetic Turbidimetric Method)	<20 EU/Sample
Bacterial Filtration Efficiency (BFE)	
Method, ASTM F2101	98.8%

TO ORDER

Re-order code

Size

S-BPC-005	5cm (2.0") Ø
S-BPC-008	8cm (3.1") Ø
S-BPC-010	10cm (3.9") Ø
S-BPC-013	13cm (5.1") Ø
S-BPC-018	18cm (7.0") Ø
S-BPC-023	23cm (9.0") Ø
S-BPC-030	30cm (11.8") Ø
S-BPC-041	41cm (16.1") Ø
S-BPC-051	51cm (20.0") Ø
S-BPC-061	61cm (24.0") Ø
S-BPC-076	76cm (29.9") Ø
S-BPC-091	91cm (35.8") Ø
S-BPC-122	122cm (48.0") Ø
S-BPC-152	152cm (59.8") Ø
S-BPC-183	183cm (72.0") Ø
S-BPC-229	229cm (90.0") Ø

Packing

S-BPC-005/S-BPC-008/S-BPC-010/S-BPC-013/S-BPC-018/S-BPC-023/S-BPC-030/S-BPC-041/S-BPC-051/:

10 pieces per sealed inner PE bag; 10 inner bags per sealed outer PE bag; Five outer bags per lined carton (500 pieces)

S-BPC-061/S-BPC-076/S-BPC-091/S-BPC-122:

10 pieces per sealed inner PE bag; 10 inner bags per sealed outer PE bag; One outer bag per lined carton (100 pieces)

S-BPC-152/S-BPC-183/S-BPC-229:

Two pieces per sealed inner PE bag; 10 inner bags per sealed outer PE bag; One outer bag per lined carton (20 pieces)

For additional information visit us at www.bioclean.com, or call us at +60 3 5569 3857/3859

Nitritex (M) Sdn.Bhd

2 Jalan Jurunilai U1/20

Seksyen U1, Hicom Glenmarie Industrial Park

40150 Shah Alam, Selangor

Malaysia

S-BPC/1809/PDS11/APAC

The Ansell logo features the word "Ansell" in a bold, blue, sans-serif font. A thick, blue, curved line underneath the text suggests a protective barrier or a seal, reinforcing the brand's association with safety and protection.

Neither this document nor any other statement made herein by or on behalf of Ansell should be construed as a warranty of merchantability or that any Ansell product is fit for a particular purpose. Ansell assumes no responsibility for the suitability or adequacy of an end user's selection of accessory for a specific application.

Ansell, ® and TM are trademarks owned by Ansell Limited or one of its affiliates. © 2012 – 2018 Ansell Limited. All Rights Reserved.